

Overview of Module 2

1. Name one protein function:

2. What is the purpose of a Western blot?

3. What is the purpose of a BCA Assay? Which metal is present in BCA reagent?

4. What is the purpose of SDS-PAGE?

5. Why do we need to transfer proteins from a gel to a membrane?

Practice Preparing a BCA Assay

1. Calculate serial dilution volumes. We will start with a 20 ug/mL BSA solution.

- a. Need 25 uL of standard for each well x 3 replicates = 75 uL per standard
- b. Always make a little extra: _____ uL per standard

BSA Concentration (ug/mL)	BSA Solution Added (uL)	Water Added (uL)	Total Volume (uL)	Final Volume (uL)
2000 ug/mL	160 uL of 2000 ug/mL	0	160	160 - 80 = 80
1000 ug/mL	80 uL of 2000 ug/mL	80	160	160 - ____ = ____
500 ug/mL	____ uL of 1000 ug/mL	____	160	160 - ____ = ____
250 ug/mL	____ uL of 500 ug/mL	____	160	160 - ____ = ____
50 ug/mL	____ uL of 250 ug/mL	____	160	160 - ____ = ____
5 ug/mL	____ uL of 50 ug/mL	____	160	160 - ____ = ____
0 ug/mL	0	160	160	160

2. Design layout for 96 well plate:

- Number of standards _____ x Number of repeats _____ = _____ wells
- Number of samples 2 x Number of repeats 2 = 4 wells
- Total number of wells = Standard wells _____ + Sample wells 4 = _____ wells
- Label the well plate below with standard and sample placement

	1	2	3	4	5	6	7	8	9	10	11	12
A												
B												
C												
D												
E												
F												
G												
H												

Copyright © 2009 Edita Aksamitiene


3. Prepare reaction buffer:

- Total number of wells _____ x 200 uL/well = _____ uL reaction buffer
- Always make a little extra: _____ uL
- Combine reagents A, B, and C in ratio 25:24:1 to make reaction buffer
 - Reagent A = (total volume of buffer _____ uL)(25/50) = _____ uL
 - Reagent B = (total volume of buffer _____ uL)(24/50) = _____ uL
 - Reagent C = (total volume of buffer _____ uL)(1/50) = _____ uL


Western Blot Hall of Shame (Challenge Question)

What do you think went wrong in the gels below?


A.

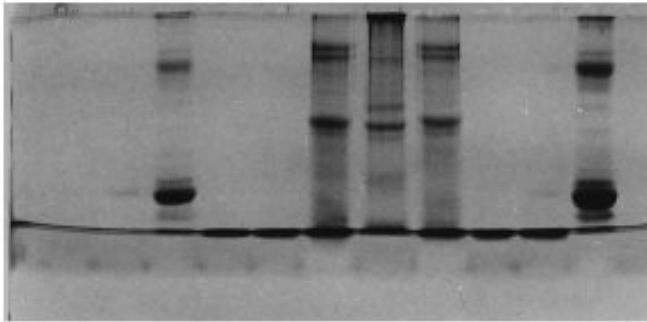


B.

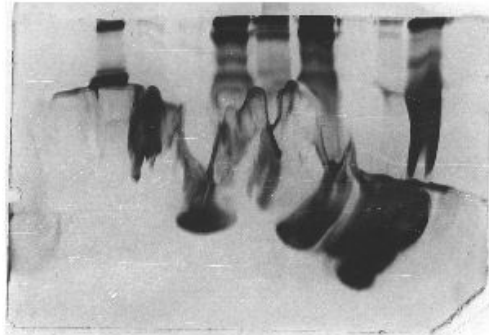


C.

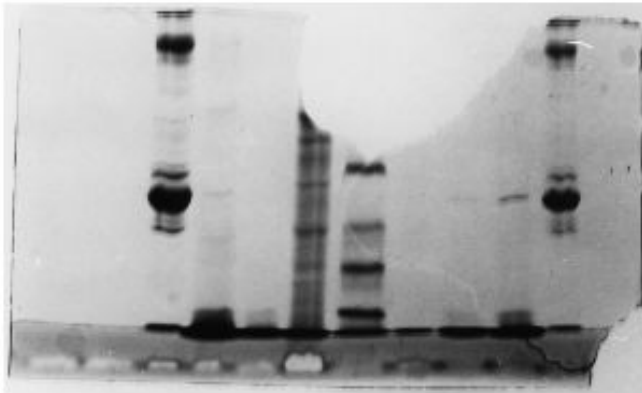


D.

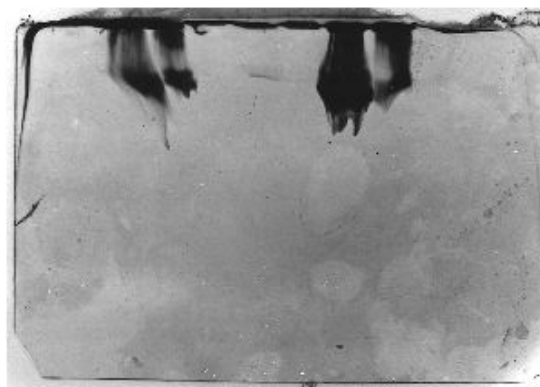

E.

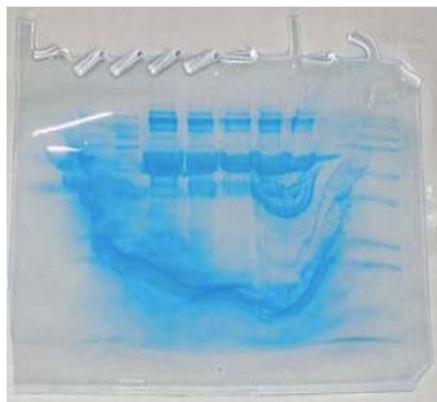

F.

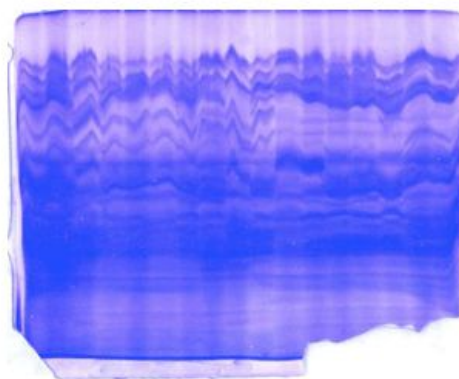

Derivation of Protein Displacement (Challenge Question)

Manipulate Newton's Second Law to demonstrate why smaller proteins travel further than larger proteins.

$$\vec{F} = m\vec{a} = m\frac{d\vec{v}}{dt} = m\frac{d^2\vec{x}}{dt^2}$$